

PIG Mercury Spill Vacuum Kit

Instructions

1. Always wear proper personal protective equipment.
2. Insert one mercury collection vial onto the nozzle of the mercury hand pump vacuum.
3. Extend flexible tubing. Place stainless steel tip into mercury spill droplets and depress the lever on the hand pump.

NOTE: For the most efficient operation, do not hold the hand pump vacuum higher than necessary above the mercury spill. Do not fill the collection tube closer than ½" below grommet. If necessary, use the scoop/spatula to help gather small mercury droplets.

4. When collection is complete, extend the tube above the pump and depress the lever on the hand pump. This will remove any remaining mercury in the tube.
5. After mercury has been collected into the mercury collection vial, carefully remove it from the nozzle of the hand pump.
6. Cut approximately 6" of butyl rubber tape and enclose grommet on mercury collection vial.
7. Place used mercury collection vial, scoop/spatula, disposable latex gloves and premoistened wipes into plastic bag provided.
8. Place hand pump into empty plastic zip closure bag provided. This will keep the potential for cross-contamination to a minimum.



DISPOSAL NOTICE: Dispose of in compliance with local, state, federal or country regulations. This product is considered nonhazardous in its purchased or unused state and requires no special disposal procedures.

SPILL KIT CONTENTS:

- (1) Mercury Hand Pump Vacuum
- (6) Collection Vials
- (1) Pair of Safety Glasses
- (6) Pair of Latex Disposable Gloves
- (6) Scoop/Spatulas
- (1) 40" Roll of Butyl Rubber Tape
- (7) Plastic Zip Closure Bags
- (6) Premoistened Wipes



New Pig

One Pork Avenue
Tipton, PA 16684-0304

newpig.com

North America: **1-800-468-4647**

Europe: **+31 (0)76 596 92 50**

China: **+86-21-400 921 5178**

PIG, PIG logo are registered trademarks in USA and other countries. See tm.newpig.com

UK: **0800 919 900**

Outside North America: **+1-814-684-0101**

165921 / 353482